

Resolution of Ether- and Diacyl-Linked Phospholipids by Differential Ion Mobility Spectrometry (DMS)

SeleXION® Technology

Cyrus Papan, Baljit K Ubhi, J. Larry Campbell and Paul RS Baker
SCIEX

The structural diversity among phospholipids is surprisingly complex, with as many as 7,000 different phospholipid molecular species in a biological sample. Often overlooked when considering structural diversity is the bond type at the *sn*-1 position of the glycerol backbone, which can be an ester bond, an ether bond or a vinyl-ether bond. Certain cell types, especially inflammatory cells and mitochondria-rich cells such as cardiomyocytes, have elevated levels of ether-linked lipids.

The Challenge:

Ether-linked and diacyl-linked phospholipids are near isobars; therefore it is necessary to resolve ether-linked and diacyl-linked phospholipid molecular species to adequately characterize their molecular species compositions. Chromatographic separation is a possibility, but due to the diversity of molecular species within a phospholipid class, this approach is not practical for profiling experiments. Even with high resolution MS, product ion analysis depends on a low-resolution isolation step that would generate a non-specific, convoluted MS/MS spectrum from the nearly isobaric compounds.

The Solution:

SeleXION technology (DMS) resolves molecules prior to MS based on their chemical properties and is effective at separating ether-linked lipids from diacyl-linked lipids in complex. Here, the presence of a carbonyl at the *sn*-1 position of the glycerol backbone changes the dipole moment and allows resolution of these phospholipid sub-classes and acquisition of non-interfered MS and MS/MS spectra for qualitative and quantitative analysis.

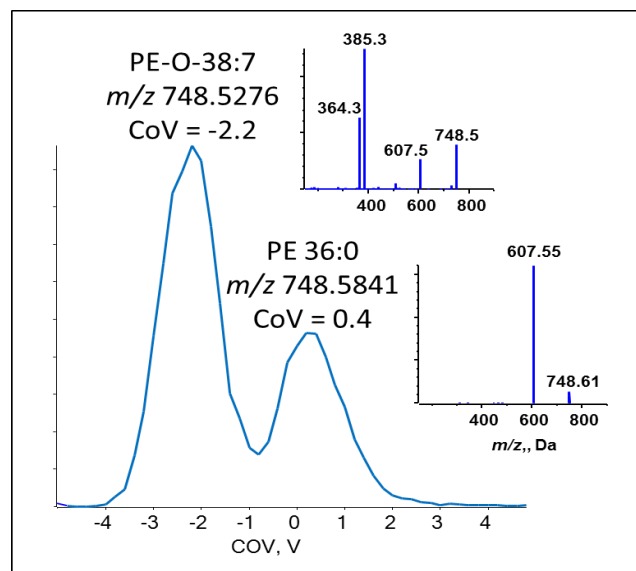


Figure 1. SeleXION Technology Resolves Ether- and Diacyl-Linked Molecular Species of Phosphatidylethanolamine. Commercially available PE extract was infused and analyzed by using a QTRAP® 6500 System equipped with SeleXION™ Technology, positive ion SIM scan mode was used to monitor m/z 748.6 while ramping the compensation voltage (CoV). Appearing in the spectrum are two distinct peaks: PE O-38:7, an ether-linked PE molecular species (i.e., PE plasmalogen with 18:1 ether-linked at the *sn*-1 position and 22:6 esterified at the *sn*-2 position of glycerol backbone), and PE 36:0, a diacyl-linked species. (Inset) Subsequent product ion analysis of m/z 748.6 using species-specific CoV values generated spectra consistent with either an ether- or diacyl-linked PE molecular species.

SeleXION Technology can resolve ether-linked from diacyl-linked PE molecular species without the need for high resolution MS or complicated chromatography.

AB Sciex is doing business as SCIEX.

© 2014 AB Sciex. For Research Use Only. Not for use in diagnostic procedures. The trademarks mentioned herein are the property of AB Sciex Pte. Ltd. or their respective owners. AB SCIEX™ is being used under license.

Document number: RUO-MKT-02-1870-A