

# Cliquid 3.4 HotFix 5 Release Notes



---

## New in Cliquid 3.4 HotFix 5

### Enhancements and Fixed Issues

Cliquid 3.4 HotFix 5 includes HotFixes 2, 3, and 4.

- This HotFix adds support for the Analyst software 1.7.4 with HotFix 1.

## Install the HotFix

| Prerequisites   |
|---|
| <ul style="list-style-type: none"><li>• The Analyst software 1.7.4 with HotFix 1 is installed.</li><li>• The Cliquid software 3.4 with HotFix 1 is installed.</li></ul> |



1. Log on to the computer as a user with Administrator privileges.
2. Stop acquisitions that are in progress, and then deactivate the hardware profile.
3. Close all applications.
4. Download **Cliquid 3.4 HotFix 5** from [sciex.com/software-downloads](https://sciex.com/software-downloads).

---

**Tip!** To prevent installation issues, save the file to the local computer, in a location other than the desktop.

---

5. After the download is complete, right-click the `Cliquid_3.4_HotFix5.zip` file.
6. Click **Extract All**, browse to and select the destination folder, and then click **Extract**.
7. Browse to and then double-click the `Cliquid_3.4_HotFix5.exe` file.
8. Follow the on-screen instructions to complete the installation.
9. Open the Cliquid software, and then activate the hardware profile. Refer to the documentation for the Cliquid software.

## Remove the HotFix

Remove the HotFix before the Cliquid software is removed.

## Cliquid 3.4 HotFix 5 Release Notes

---

**Note:** For SCIEX 4500MD systems and Citrine systems with ICB-5, this HotFix should not be removed because ICB-5 is supported only if this HotFix is installed.

---

1. Log on to the computer as a user with Administrator privileges.
2. In the Analyst software, deactivate the hardware profile.
3. Close the Cliquid software.
4. Close all applications.
5. Open Control Panel, and then open Programs and Features.
6. Select **Cliquid 3.4 HotFix 5**, and then click **Uninstall**.
7. Obey the on-screen instructions.

The HotFix is removed from the program list, and any files that were changed by the HotFix are reverted to the Cliquid 3.4 with HotFix 1 version.

---

**Note:** Any HotFix that was removed during installation of the Cliquid 3.4 HotFix 5 is not restored. It must be installed again after HotFix 5 is removed. A compatible version of the Analyst software must be installed to make sure that the Cliquid software 3.4 operates correctly.

---

8. Close Control Panel.
9. Remove the Analyst software 1.7.4 with HotFix 1, and then install a version of the Analyst software that is compatible with the Cliquid software 3.4 with HotFix 1.
10. Open the Cliquid software, and then activate the hardware profile. Refer to the documentation for the Cliquid software.

## Updated Files

The Cliquid 3.4 HotFix 5 updates the following files:

**C:\Program Files (x86)\AB SCIEX\Cliquid\Client**

- Cliquid.exe
- Cliquid.Resources.dll

**C:\Program Files (x86)\AB SCIEX\Cliquid\Server**

- Analyst.reg
- tech.reg

## Known Issues and Limitations

- N/A

## Contact Us

### Addresses

|                               |   |
|-------------------------------|---|
| <b>SCIEX<br/>Headquarters</b> | AB Sciex LLC<br>250 Forest Street<br>Marlborough, MA 01752<br>USA |
|-------------------------------|---|

## Customer Training

- Global: [sciex.com/contact-us](https://sciex.com/contact-us)

## Online Learning Center

- [SCIEX Now Learning Hub](#)

## SCIEX Support

SCIEX and its representatives have a global staff of fully-trained service and technical specialists. They can supply answers to questions about the system or any technical issues that might occur. For more information, go to the SCIEX website at [sciex.com](https://sciex.com) or use one of the following links to contact us.

- [sciex.com/contact-us](https://sciex.com/contact-us)
- [sciex.com/request-support](https://sciex.com/request-support)

## Cybersecurity

For the latest guidance on cybersecurity for SCIEX products, visit [sciex.com/productsecurity](https://sciex.com/productsecurity).

## Documentation

This version of the document supersedes all of the previous versions of this document.

To see this document electronically, Adobe Acrobat Reader is required. To download the latest version, go to <https://get.adobe.com/reader>.

To find software product documentation, refer to the release notes or software installation guide that comes with the software.

## Cliquid 3.4 HotFix 5 Release Notes

---

The latest versions of the documentation are available on the SCIEX website, at [sciex.com/customer-documents](https://sciex.com/customer-documents).

---

**Note:** To request a free, printed version of this document, contact [sciex.com/contact-us](https://sciex.com/contact-us).

---

---

This document is provided to customers who have purchased SCIEX equipment to use in the operation of such SCIEX equipment. This document is copyright protected and any reproduction of this document or any part of this document is strictly prohibited, except as SCIEX may authorize in writing.

Software that may be described in this document is furnished under a license agreement. It is against the law to copy, modify, or distribute the software on any medium, except as specifically allowed in the license agreement. Furthermore, the license agreement may prohibit the software from being disassembled, reverse engineered, or decompiled for any purpose. Warranties are as stated therein.

Portions of this document may make reference to other manufacturers and/or their products, which may contain parts whose names are registered as trademarks and/or function as trademarks of their respective owners. Any such use is intended only to designate such products as those manufacturers' products and does not imply any right and/or license to use or permit others to use such manufacturers' and/or their product names as trademarks.

SCIEX warranties are limited to those express warranties provided at the time of sale or license of its products and are the sole and exclusive representations, warranties, and obligations of SCIEX. SCIEX makes no other warranty of any kind whatsoever, expressed or implied, including without limitation, warranties of merchantability or fitness for a particular purpose, whether arising from a statute or otherwise in law or from a course of dealing or usage of trade, all of which are expressly disclaimed, and assumes no responsibility or contingent liability, including indirect or consequential damages, for any use by the purchaser or for any adverse circumstances arising therefrom.

(GEN-IDV-09-10816-E)

For Research Use Only. Not for use in Diagnostic Procedures.

Trademarks and/or registered trademarks mentioned herein, including associated logos, are the property of AB Sciex Pte. Ltd., or their respective owners, in the United States and/or certain other countries (see [sciex.com/trademarks](https://sciex.com/trademarks)).

AB Sciex™ is being used under license.

© 2025 DH Tech. Dev. Pte. Ltd.

